Golden Rules For CHEMICAL FUME HOODS

- 1. Make sure the hood is working.
- 2. Keep the sash down except when working in the fume hood.
- 3. Raise the sash only the minimum needed to perform the work.
- 4. Work six to eight inches inside the plane of the sash.
- Keep hood area clear and rear vent unobstructed.
- 6. Raise equipment off hood deck. (Use PVC pipe donuts, not wood.)



- 7. Remove flammables and unnecessary stuff.
- 8. Understand the limitations.
- 9. Check / inspect operation at least annually, sash cables every 2-3 years.
- 10. Wear proper personal protective equipment (PPE).
 - cables and

11. Run cords, cables and tubing under the air foil.

12. Move arms slowly in and out of the hood.



More at **LabSafety.org.** The Laboratory Safety Institute is a non-profit organization committed to making health, safety and the environment an integral and important part of education, work, and life.



I

